

FastGene® CapEasy



Check it out!

The FastGene® CapEasy (FG-CDC01) was created to ease up your daily labwork. Capping or decapping the lid of an 8- and 12-well strip is complicated and often leads to the rupture of the lid. This can lead to a cross-contamination between the wells making a repetition of the whole experiment necessary.



Cat.No.: FG-CDC01

PCR-Tube recommendation

- FG-028 combined with:
 - FG-008FC (flat caps)
 - FG-008DC (domed caps)
- FG-016FC (flat caps)
- FG-016DC (domed caps)

This can be avoided by using the FastGene® CapEasy. The uniformly distributed pressure ensures the sealing of all wells and it doesn't matter if it is a domed lid or a flat lid. The handling of the FastGene® CapEasy is, as the name already says, very easy:

Procedure

Close all the lids at once!

Place the strip in the green rack. Fix the position of the caps lightly on top. Close the lid without using too much strength.



Close the lid easily!



Open all the lids at once!

Place the strip with the caps in the green rack. Position the lid lightly on the caps. Slide the comb to the tubes until a click-sound is heard.

